SUPACARB CP

High Performance Carbon Block

SupaCarb CP offers the very best in activated carbon technology. The cartridge is manufactured using a unique sintered carbon that is specifically designed for maximum decolourisation and organic contaminant reduction. It offers far superior adsorption levels when compared to other carbon block cartridges.

SupaCarb CP is an extremely economical and effective solution to many common liquid treatment applications where the reduction of colour, organics and other impurities is required. The carbon block is made using a high purity micro-porous, acid washed, steam activated carbon and is particularly suited to pharmaceutical applications.



- Unique sintered block using both granular and powdered carbon
- Spun bonded polypropylene prefiltration layer
- Available from 5mm segments to 1018mm (40") long cartridges, suitable for laboratory, pilot plant and production scale operations.



The SupaCarb CP carbon block uses a unique sintering process that combines both granular and powdered activated carbon. This results in the highest available surface area for adsorption of contaminants. The carbon block is protected by a spun-bonded, pure polypropylene fibre prefiltration layer, to produce a unique high performance module, offering superior on-stream life.

SupaCarb CP cartridges are manufactured using the latest thermal welding technology, eliminating bypass commonly associated with adhesive bonded carbon modules. This results in a robust and durable filter assembly. They are suitable for a wide range of process conditions, including use in the harsh operating conditions encountered in pharmaceutical API plants.

Features and Benefits

- Combination of both granular and powdered activated carbon offers superior contaminant adsorption levels
- Low volume pre-use flush required compared to lenticular cartridges
- · High active carbon content per module results in the use of smaller installations, less space required and lower costs
- Convenient low-weight cartridge format is easy to handle and modules can be changed-out without using lifting gear, special tools or other apparatus
- High purity micro-porous steam activated powdered carbon is particularly suited to removal of coloured impurities from pharmaceutical product and solvent streams
- Graded density prefiltration layer protects carbon block offering extended life
- Thermally bonded construction eliminates fluidising, channelling and bypass
- All polymeric support structure offers excellent chemical compatibility
- Each cartridge is embossed with an identification number and fully traceable

Industries and Applications

Pharmaceutical and Fine Chemicals

Food and Beverage

Water Treatment

Chemical Manufacturing

- Removal of coloured impurities from product and solvent streams
- Colour and odour reduction in beverages
 - Removal of chlorine, ozone, odour, colour and trace organic compounds
- Removal of catalysts and wastewater treatment



SupaCarb CP Technical Data

Dimensions

Outside Diameter: 66.5mm Inside Diameter: 27mm

Sterilisation and Sanitisation*1

Steam: 121°C for 15 mins (20 cycles) Hot Water: 90°C for 30 mins (0.2 bar Δp max) "Applies to single open end cartridges only. For all steaming and hot water applications, the Reinforced Polypropylene End Cap option must be used .

Maximum Operating Conditions

80°C Temperature:

Recommended change-out

differential pressure: 2.5 Bar

Recommended flow rate: 2.0l/minute per 10" filter

Product validation guide available on request. All SupaCarb CP cartridges are manufactured under strict control with batch number identification, giving full traceability on all components.

Ordering Guide

| 06CP | 005 - | 20 | 3 | E | Α | |
|------------------------------|---------------|---|---|--|------------|--|
| Carbon/Media | Micron Rating | Length | Connections | Seal | Branding | Options |
| 06CP - Carbon Sintered Block | 005 - 5.0μm | 09 - 249mm 20 - 510 30 - 764 40 - 1018 | 0 - DOE A - Code A B - Code B*2 S - Code S 2 - Code 2 3 - Code 3 7 - Code 7 8 - Code 8 | P - PE Gasket B - Buna E - EPDM F - FEP/Silicone (SOE Only) S - Silicone V - Viton | A - Amazon | G - Reinforced Polypropylene End Cap |

Example: $06CP005-203EA = Carbon sintered block media, 5.0 \mu m rating, double length 20" long, code 3 connections with EPDM seal. *2 Code B - to fit Amazon 50 Series housings only.$

Amazon Filters Ltd.

Albany Estate, Camberley, Surrey, GU16 7PG, ENGLAND

Tel: +44 (0) 1276 670 600 Email: sales@amazonfilters.co.uk Web: www.amazonfilters.com